



Shortcuts of JENOPTIK GRYPHAX® software

JENOPTIK GRYPHAX® software is workflow-optimized and very easy to use. Therefore, JENOPTIK GRYPHAX software includes a lot of time-saving keyboard shortcuts.

We all want to boost our productivity, so we recommend printing out or saving this list of keyboard shortcuts to enhance your daily work!

Keyboard button:	Shortcuts for:
F1	Open help
F2 ENTER	Start capturing in "Capture Mode" (single shot, time-lapse, video)
F3	Return to " Live " mode
F4*	Open Individual Save dialog
F5	Activates deactivates live image freeze
F11	Switch between " Presentation view " and "Standard view"
DEL	Delete selected images from "Gallery" or marked drawings (measurement annotation)
Z	Switch zoom between "1:1" mode and "Fit to screen" mode
CTRL** + P	Print selected image(s) from "Gallery"
CTRL** + A	Select all (all thumbnails from "Gallery")
CTRL** + G	Open close " Gallery "
CTRL** + O	Open " Storage Options " to change destination folders for gallery (images, videos)
CTRL** + T	Open close " Toolbar "
CTRL** + D***	Open close " Treeview " to change and manage media destination folder fast & easy
G	Open close " Grid " view
ESC	Leave "Presentation view" remove drawings during creation (measurement annotation)
Arrow up down	Gallery thumbnail up down
Page up down	Gallery page up down
Home	Jump to "Gallery" first page
End	Jump to "Gallery" last page
SPACE BAR	Activates " Magnifier " or " Hand tool "
+ -	to zoom IN or OUT while "Magnifier" is active

*valid as of GRYPHAX version 2.1.0 | **for Mac users: CMD instead | ***valid as of GRYPHAX version 1.1.6



MORE LIGHT

Learn more about JENOPTIK GRYPHAX® software and the easy workflows and tools.

Watch our video tutorials.

[VIDEO TUTORIALS](#)

HELPFUL? We appreciate your feedback.

[FEEDBACK](#)

We looking forward to hearing from you and remain sincerely,

Your JENOPTIK GRYPHAX® Team

GRYPHAX® SERIES

Release date: 18/06/2019

JENOPTIK | Light & Optics
JENOPTIK Optical Systems GmbH
Goeschwitzer Strasse 25
07745 Jena, Germany